

<b>Title:</b>	<b>Saving Permanent Changes of Configuration DMs</b>	<b>ID:</b>	
		0178	
<b>Date in:</b>	<b>Response:</b>	<b>Model:</b>	<b>Author:</b>
2003-09-04	2003-09-04	-	CMa

## Q:

Is there a way to permanently set the number of hotel stacks, shelves/hotel, and shelf pitch (DM29, DM25, DM23) permanently in the firmware like we do with the z-travel, turn axis, etc. variables that we save with the "Thermo Updater" program?

Can you provide us with a tool to do this?

Also, for the open air carousel, what step size should we use for the deep well plate (DM23)?

## A:

There are two proposals:

1. You send us the serial number of the system ("RD DM109") and the constants to put in and we will provide you a COD-File and a download program.
2. Our ActiveX library supports a configuration feature that allows you to edit the DM-settings. A call of this function will update the DM's

The open carousel has 788 vor 23mm -> for DWP there would be 50mm = 1713 steps.