

Title:	What is second d-sub connector on the back			ID:
				0254
Date in:	Response:	Model:	Author:	
2005-02-08	2005-02-08	-	C <i>Ma</i>	

Q

Thanks a lot by your information. I'm sure it will be useful for us. I have only one additional question:

The optional port that appears beside the RS232 controller like on the Cytomat incubators, to give information about the Temperature, RH% and CO2.

A

The second port on Evotec Systems is the CAN through on other units its the BCR port.

On the cytomat you had two ports because on these units we retro-fitted the handling in an standard incubator. Liconic combined climate and handling control. So you may request the information through the same port.

To read

- Temp. RD DM982 (actual) RD DM890 (setting)
- RH RD DM983 (act.) RD DM893 (set.)