

Title:	CO2 not controlling to set-point	ID:	
		0260	
Date in:	Response:	Model:	Author:
2005-03-08	2005-03-08	-	WOp

CO2 not controlling to set-point

1. check if the set-point value is correctly programmed (RD DM894)
2. check if the correct gas is connected to the unit
3. check the incoming CO2 pressure
4. check if the sensor pump is ON (RD 701) PLC channel 701
5. check if solenoid sensor is turning On and Off (RD 700) PLC channel 700
6. check delay timer (T92) if low set to (WS T92 65000)
7. reset T92 if needed (RS)