

Title:	Error 18	ID:	
			0347
Date in:	Response:	Model:	Author:
2009-12-08	2009-12-08	STX220	CMa

Q:

STX220 returned 10258 to DM200. ErrConv software says it is Err18.
Remote Operation V0710 doesn't have the error code.

Please advise what error it is.

A:

Error 18 is the carrousel positioning error. Error 18 was implemented in V7.1.16. By mistake Error 18 was also assigned to Shovel Retract Timeout in V7.1.17. The bug has been fixed in V7.1.25. where Shovel Retract error has become Error 20