

Title:	Difference between SA and BT	ID:	
		0080	
Date in:	Response:	Model:	Author:
2002-03-09	2002-03-09	-	CMa

Q:

What is the difference between StoreX-..SA and StoreX-..BT?

A:


SA stands for StandAlone while BT stands for BenchTop .

The StandAlone stands on the ground and has the automatic Gate at ~1000mm above ground (High Gate). The Gate opens downwards. The BenchTop is placed on a table and has its automatic Gate ~200mm above the table surface (Low Gate). The Gate opens upwards. We recommend to use StandAlone units whenever possible since:

- StandAlones creates new table space rather than use up table space
- StandAlones are easier to place and / or move (use fork lift or a trolley)
- StandAlones are easier to upgrade from a smaller to a larger unit
- StandAlones allow free access from the top since the Gate moves downwards

BenchTops are mainly used when there is an existing table.

Example: STX40-..SA

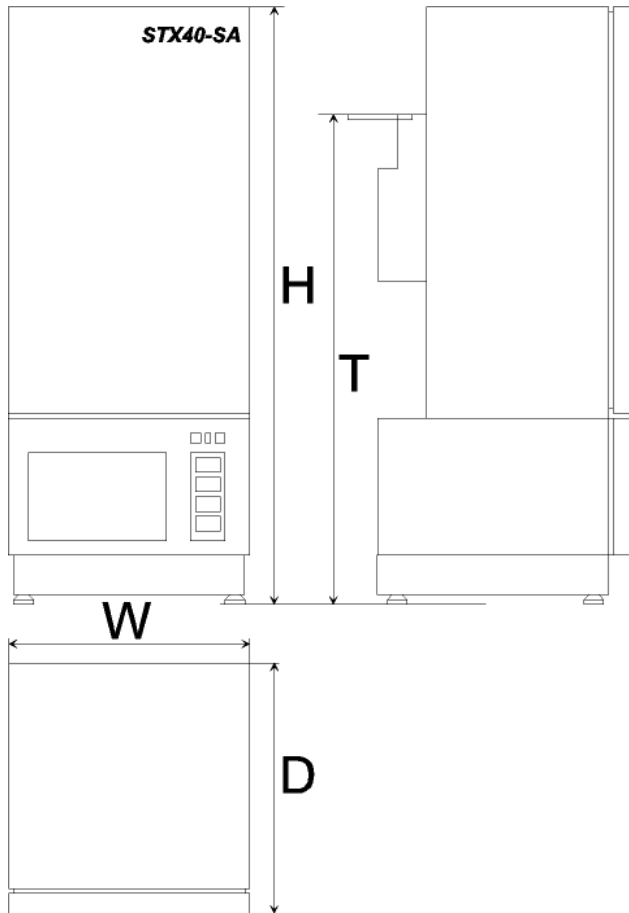

Example: STX40-..BT

