

|                 |                         |               |                |
|-----------------|-------------------------|---------------|----------------|
| <b>Title:</b>   | <b>DWP MTP Handling</b> | <b>ID:</b>    |                |
|                 |                         | 0095          |                |
| <b>Date in:</b> | <b>Response:</b>        | <b>Model:</b> | <b>Author:</b> |
| 2003-05-24      | 2003-05-24              | -             | CMa            |

**Q:**

I would like to confirm that all incubator mechanisms are able to handle both deepwell plates and microplates by just changing the racks, pitch and transfer station position.

**A:**

Unfortunately they are not. There is a mod-kit that increases the height of the shovel. Ideally this should be specified at the time of order.

The new handling used in the STX500, STX1000 and STX220 can handle both DWP and MTP. The plan is to introduce this new handling to the STX40 series this year.