

Title:	Error Code	ID:	
			0098
Date in:	Response:	Model:	Author:
2003-03-05	2003-03-05	-	CMa

Q:

I got an error code that's not in the manual.
I issued RD 1814 and I got back 1 indicating that there's an error.
So I issued RD DM200 and I got back 08961.
What is 08961 mean? It's not in any of the documents that I have.
Could you please explain this.

A:

8961 is Hex \$2307. Obviously you did an Unload (2) and your handler got blocked when extracting the shovel at the transfer station (3) and the Gate did not open.. I added the command set. Your questions concerning the ActiveX will be answered by Maxim.