



<b>Title:</b>	<b>Cooling Ventilator Motor Connection</b>	<b>ID:</b>	
		0054	
<b>Date in:</b>	<b>Response:</b>	<b>Model:</b>	<b>Author:</b>
2007-11-17	2007-11-17	STP1000	CMa

## Connection Terminals Cooling Ventilator STP1000 / STX1000-DF

There are two types of motors used for cooling in the STX1000-DF. Systems built prior 2007 use the Karl-Klein motor (black, aluminum).

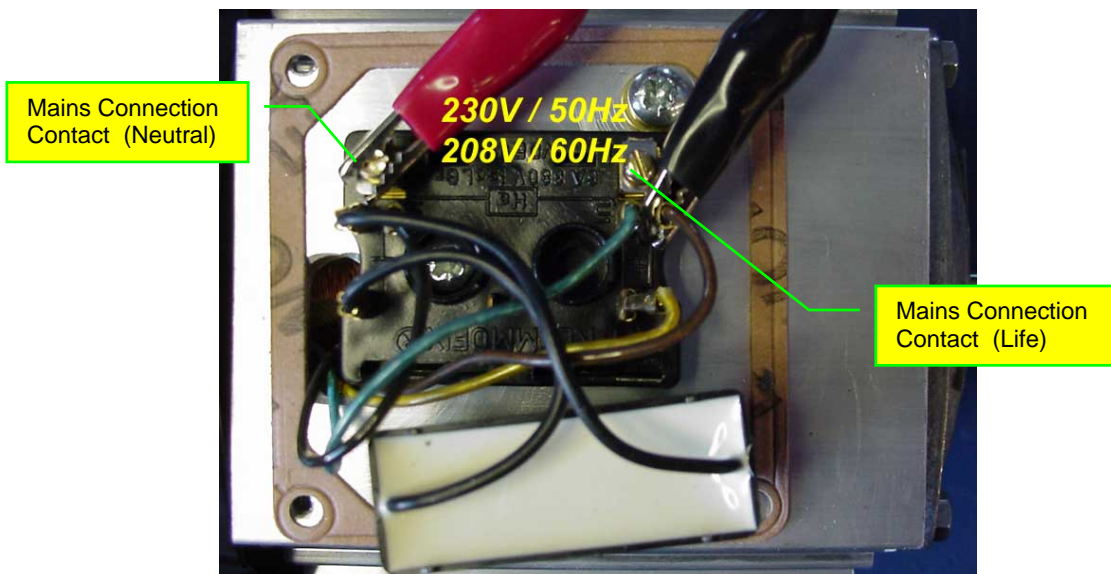

Systems built 2007, Dec. 2006 and later use the new style motor (blue) that has less heat dissipation.

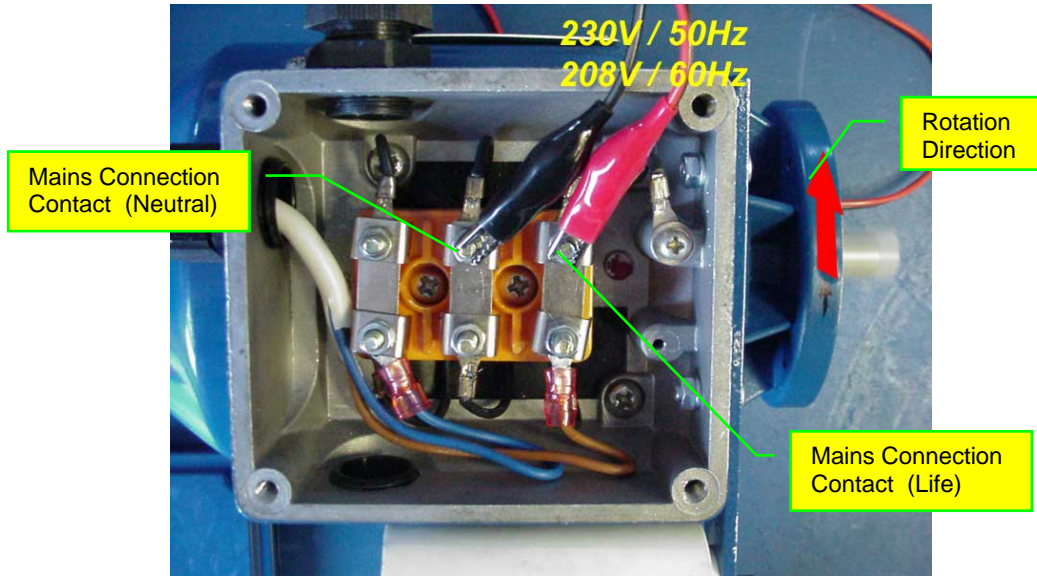

Connecting Life to the left most terminal will reverse rotation direction.